


CASY[®]-Technology


Fermentation Monitoring Bacteria (Escherichia coli)

Control of batch fermentation starting with a stationary overnight culture.


a) Shift in size distribution, exponential / stationary

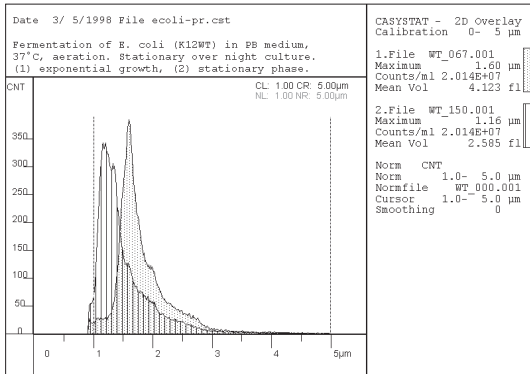

b) Change in mean cell volume during the run

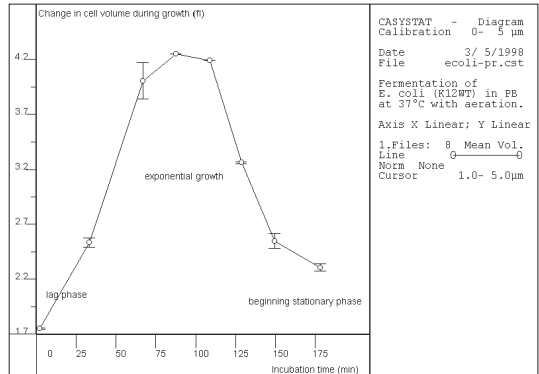

c) Proliferation

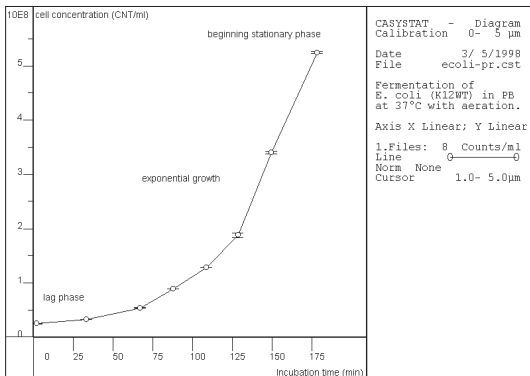

d) Analysis of the generations generated per time

Generation Times

Printing On

Gen. Time	# Generations
Set 1: 3.406e+01	2.202e+00
Set 2:	
Set 3:	
Set 4:	
Set 5:	

x-Range

From	To
100.00	175.00

By CASY[®] Model TTC using a 45 µm capillary

Copyright[®] by innovatis AG